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-turn the button on

- Open the machine cover by pressing on the door gently and direct the locker toward opening.
- Adjust the water level of the machine using distilled water.
- Put plastic or glass material to be autoclaved after put autoclave taps on each of them.
- Close the machine by direct the locker toward closing.
- Connect the electricity and press on start button .

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تعليمات انتهاء عمل الجهاز :-
- When autoclaving completed and temperature reach 80oc press on stop button and press gently on the door and direct the locker toward opening

اجراءات السلامة والامان :-
- Do not run the machine if there is problem in any parts.
- Open the autoclave slowly
- Wait until all compressed air completely free.

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- Open the machine cover by moving 6 round keys anti'clockwise.
- Adjust the water level of the machine using distilled water.
- Put plastic or glass material to be autoclaved after put autoclave taps on each of them.
- Close the machine by moving 6 round keys clockwise.
- Connect the electricity, when temperature will be 121c and pressure will be 0.1 m pa the valve will turned on automatically then adjusted to set time(20min for plastics,30 min for glasses by timer)

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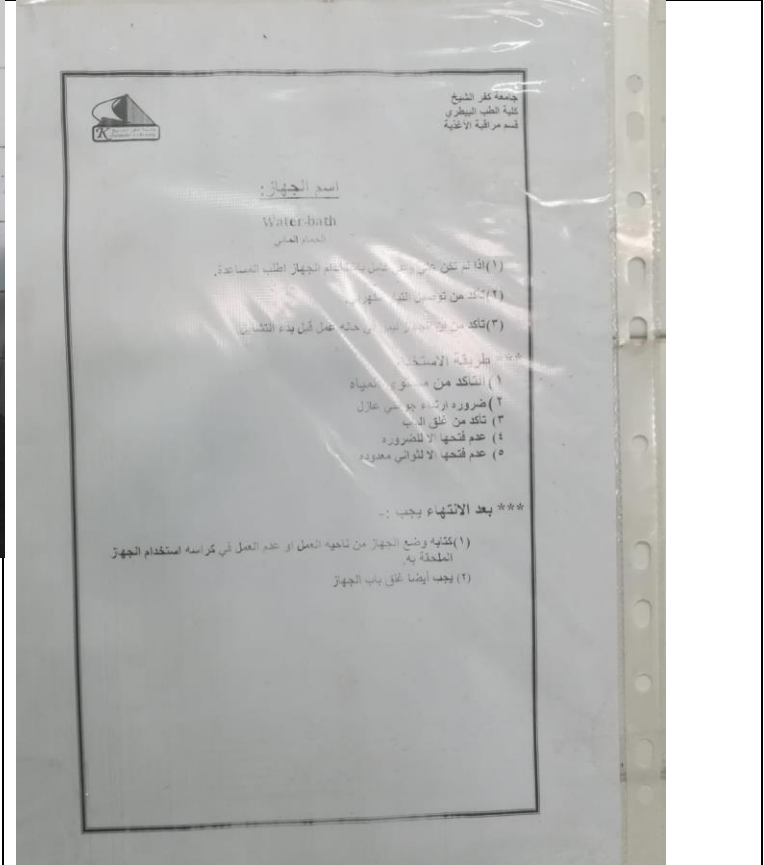
تعليمات إنهاء عمل الجهاز :-
- Turn off the machine ,wait until all air out of through the valve, observe the arrow on the machine counter till pressure will be zero open the cover and take the material, observe the autoclave taps turned back or not as a marker for successful autoclaving ,,put the cover on its proper place

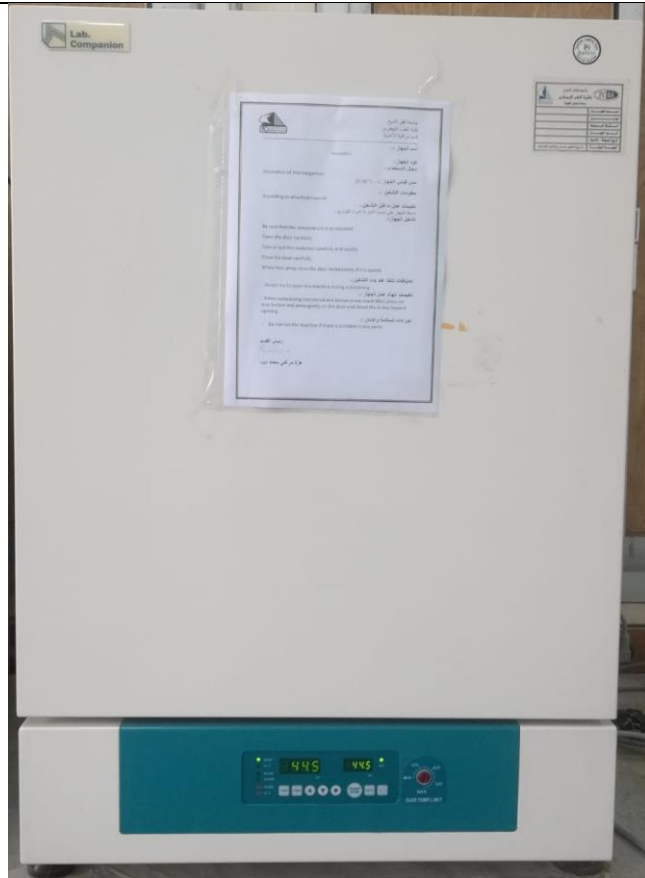
اجراءات السلامة والامان :-
- Do not run the machine if there is problem in any parts.
- Open the autoclave slowly
- Wait until all compressed air completely free.

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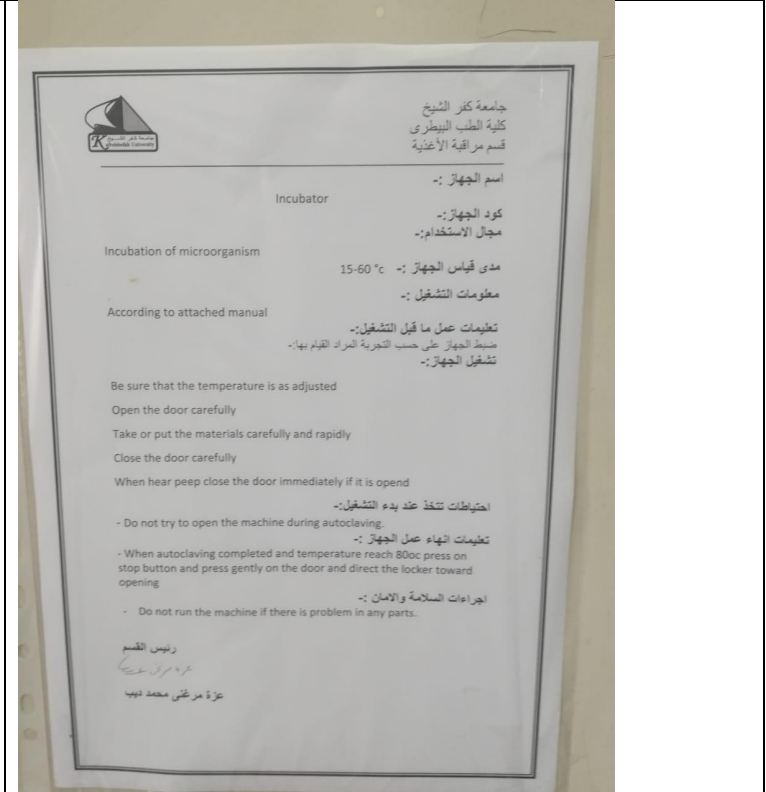
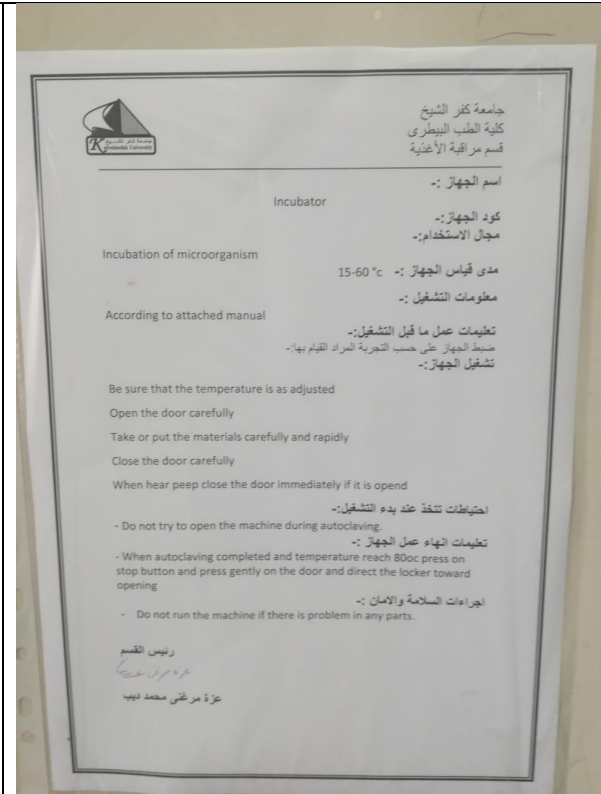
Be sure that the temperature is as adjusted
Open the door carefully
Take or put the materials carefully and rapidly
Close the door carefully
When hear peep close the door immediately if it is open

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Turn the key to 0 position
Open the machine door on the upper part by the two keys on it.
Clean the internal surface of machine using STAVELON
Turn the light on by light key

احتياطات تتخذ عند بدء التشغيل :-
Remove the working response
Clean the surface
Turn the key to position 0
Close the door of machine
turn the uv on 10 min .
Remove the electricity plug .

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- Carefully read and follow the instructions.
- Pay special attention of the load.
- Trial runs and operation under DIN protection class 0 only with continuous observation and monitoring.
- Materials sensitive to temp and valuable items must only be placed in oven

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- Connect electricity
- Open the door of the oven and insert the instrument needed to be -- sterilized.
- Close the door.
- Adjust the temperature.
- Adjust time.

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- The operation is continually observed and monitored by an operator.
- The oven is fitted with over temp protection to DIN protection class2 (adjust templimiter) class3.1(adjustable temp monitor)

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- Do not run the machine if there is problem in any parts.
- Open the oven slowly

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